

PRODUCT INFORMATION

MycoDecon

Art.-Nr. 34206

Product Description:

General Highly effective disinfectant, particularly active against mycoplasmas, but also against bacteria, virus and fungi.

Features

- Ready-to-use
- Alcohol based
- Non-corrosive
- Non-carcinogenic

Storage Store at + 15 °C to + 30 °C.

Application MycoDecon is special spray formulation for the use on any laboratory surface and equipment, including benches, incubators, storage boxes of cells, containers for liquid nitrogen.

Decontamination of surfaces:

- Spray MycoDecon directly on the contaminated surface until it is moistened completely.
- After 5 min dry the surface with a clean paper towel.
- **IMPORTANT:** Remove completely, do not leave any residue.

Please note: May cause discoloration of sensitive surfaces. Test on small part of the surface prior to use.

Decontamination of laboratory devices:

- Apply MycoDecon on a paper towel and clean all surfaces of the devices thoroughly.
- Dry with a clean paper towel.

Version 07/24