

PRODUCT INFORMATION

SERVA Unstained Protein Standard 6.5 - 97 kDa

Cat. No.: 39214


PRODUCT DESCRIPTION

SERVA Unstained Protein Standard is a ready-to-use protein marker that contains 7 proteins ranging from 6.5 to 97.4 kDa.

Application SDS-PAGE

Features

- Number of bands: 7
 Phosphorylase B (97.4 kDa)
 Albumin bovine (BSA, 67 kDa)
 Ovalbumin (45 kDa)
 Carbonic anhydrase (29 kDa)
 Trypsin inhibitor, soybean (21 kDa)
 Cytochrom C (12.5 kDa)
 Trypsin inhibitor, bovine (6.5 kDa)
- Ready to use solution
- Concentration:
 1 mg/ml in Tris/Glycine SDS sample buffer

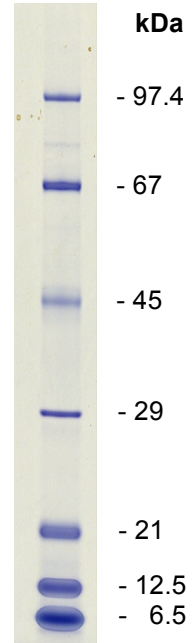

Storage

-20 °C, please aliquot in a small volume for prevention of repeated freezing & thawing.

Usage information

- Recommended loading volume for mini gel (10 x 8 cm, 0.75 or 1 mm thick)
- Coomassie staining: 5 µl
 - Silver staining: Dilute 1:5 in 1x Laemmli sample buffer and load 5 µl