

## PRODUCT INFORMATION

**Grotanol® FF 1 N**

**Cat. No. 43440**

---

### Product description:

**General**

Microbiocidal system cleaner for cleansing of circulation systems.  
Sufficient for 5L cleaning solution.

---

### Use:

- **Preparation:**  
Dilute 3% Grotanol in distilled water. Fill chiller to the maximum mark.
  - **Cleaning Procedure:**  
Turn off temperature control of the chiller or set to room temperature. Start chiller and let circulate for 24h. Afterwards, remove the solution and fill the system with distilled water. Let circulate for 5 minutes, discard the water and repeat until it is free of any impurity.
  - **Control:**  
Proof flow rate and repeat the process if the rate is still low.
- 

### Storage conditions:

- +15 °C - +30 °C
- 

Ver 0423