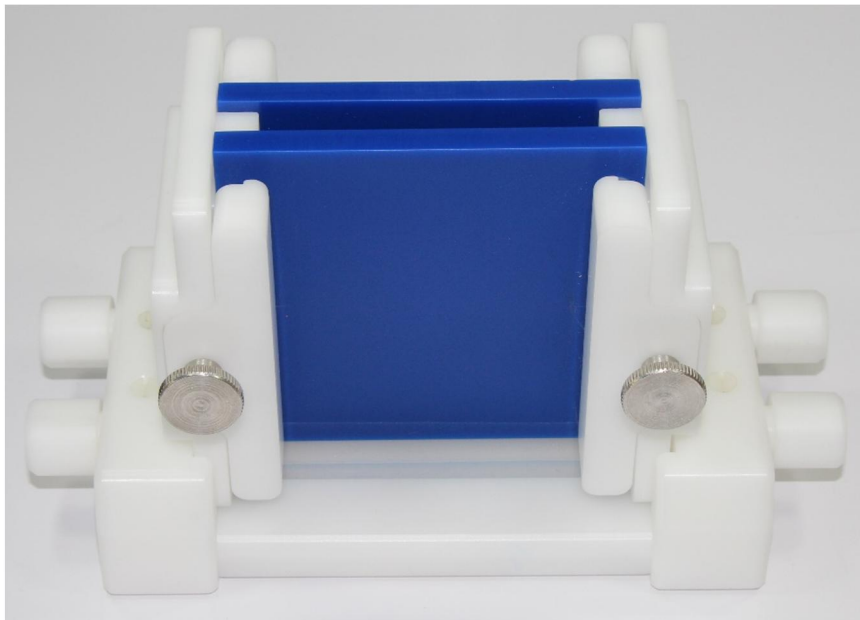
# BlueVertical™PRiME™

Gelgießstand, Casting Stand

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## General information

Keep the instruction manual close at hand nearby your device. We structured this document in a way that you can refer to the desired information via the index. We recommend to read all the chapters to obtain detailed explanations and operations notes. Aim of this instruction manual is to make the operation accessible to you with a language easy to understand. Get in touch with our technical service should any questions arise or explanations being unclear (contact details, see rear page).

By using the packing list, please check after unpacking, if all parts of the device are complete and the device is undamaged. If this is the case, please inform SERVA Electrophoresis GmbH at once.

**The warranty period is 12 month and starts at delivery. We ask for setting the packing material aside until warranty period is expired.**

## Our target group

The SERVA BlueLine devices are intended for laboratory use. English is considered as the global language of science. The target group of this instruction manual is skilled laboratory staff. Among them count workers-in-training (e.g. trainee, apprentice, diplomate and postgraduate) after briefing as well. Local safety regulations (e.g. biological, chemical, radioactive and medical risk- und security level) cannot be considered in this manual. The liability in these issues lies with the user.

## SERVA assistance

Assistance	Contact
Sales Team (Germany)	Contact your local sales representative to receive product information, to arrange demonstrations or to inquire quotations and product samples.
Distributor (International)	Outside Germany, please contact your local distributor for product information, pricing and inquiries. The contact details of our worldwide distributors are listed on our website.
Customer Care	Our sales department is informing you about prices, inquiries and shipment. The contact details are listed on the rear page of the manual.
Technical Service	Our Technical Service is your contact point for technical and scientific questions about our products. The contact details are listed on the rear page of the manual.
Product Specialist	Our Product Specialists for individual applications are glad to assist you with improving your method and solve problems. Ask our Technical Service for the contact.
Download Center	In our Download Center you find all Manuals, ApplicationNotes and TechNotes regarding our products. Downloadable are brochures, the online catalog and various certificates as well. Additionally on the single product pages in the online shop, MSDS and the specifications are deposited.
Webinar Archives	Inform yourself about our future webinar dates under Events on our website. Recorded videos of previous webinars can be downloaded from our Webinar Archives.

## General lab safety information

- In the lab, wear gloves, lab coat and safety goggles. Do **not** wear jewelry or watches.
- Avoid contact with mouth, nose and eyes by hand before having your hands washed. Treat and dress small wounds sufficiently.
- Before leaving the lab, take your safety equipment off and wash your hands thoroughly with soap.
- Change the gloves frequently and remove them before using a telephone, a light switch or a writing utensil.
- Clean your equipment, lab bench and devices frequently and directly after contamination with a mild soap and disinfectant.

## Hazards

If the user follows the safety regulations as described below, the SERVA BlueLine devices are designed in a way that safe working is ensured. Any warranty claim will be voided if the device or its spares are changed and modified by an unauthorized person. Working contrary to regulations will void warranty claim additionally.

## Cleaning and disinfection

Clean the SERVA BlueLine devices before first start-up and then frequently with a mild detergent (0.1% SDS solution) followed by distilled water as follows:

First, disconnect the device from the grid before start cleaning. Use lintless cloth soaked in water only. To remove tenacious dirt, lintless cloth soaked in 0.1% SDS solution can be used before. Never use aggressive cleaner or solvents. Let everything dry before reconnecting and start working. Parts of the device that were in contact with other liquids than water have to be cleaned after each usage to avoid salting-out and encrusting.

If contaminated, choose the disinfection method fitting your local regulations and guidelines. First, disconnect the device from the grid and let the device cool down before start disinfecting. After disinfection clean as described before.

If the device needs to be shipped back to SERVA, perform cleaning and disinfect if necessary. Document this on the Decontamination Certificate (Download Center on [www.serva.de](http://www.serva.de)) and enclose it in the shipping box.

## CE Certification

**Important:** This SERVA product is designed and certified to comply the safety guideline EN61010-1.

CE certified products are safe in use if utilized as described in the manual. The device has not to be modified. After any modification, warranty and CE certification becomes null and void. Additionally, these modifications may represent a potential source of danger.

SERVA is not responsible for any harm or injury that was caused by device modification, improper use or unspecified applications.

## Imprint

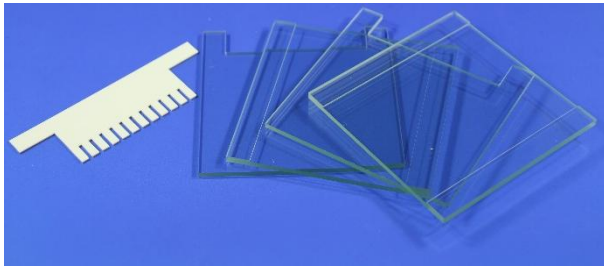
Parts of this manual neither may be changed nor used in any other form without permission in written form on the part of SERVA Electrophoresis GmbH. We reserve the right to modify our products and documents anytime. We assume no responsibility for errors, damage and injuries caused by unspecified applications or noncompliance with the advised safety regulations.

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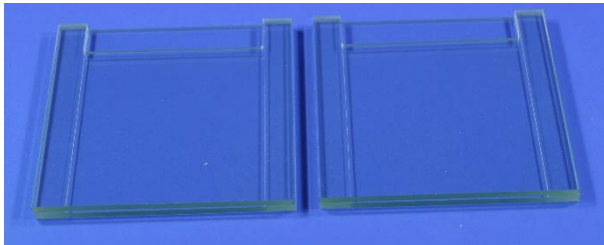
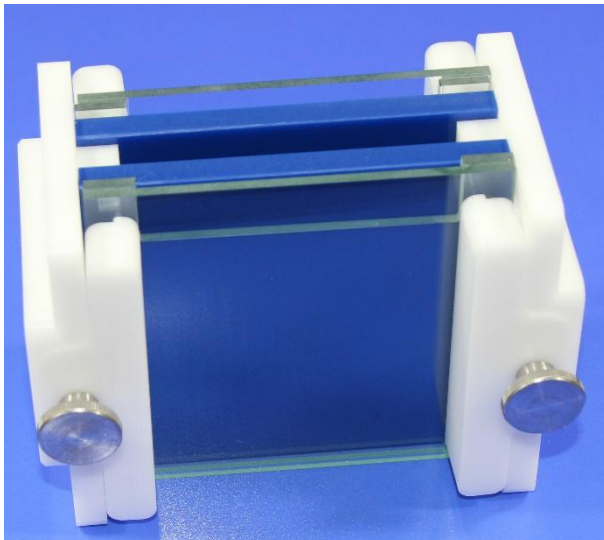
Place the plain plates with the spacers towards you onto the bench. Superpose the notches with the spacers.

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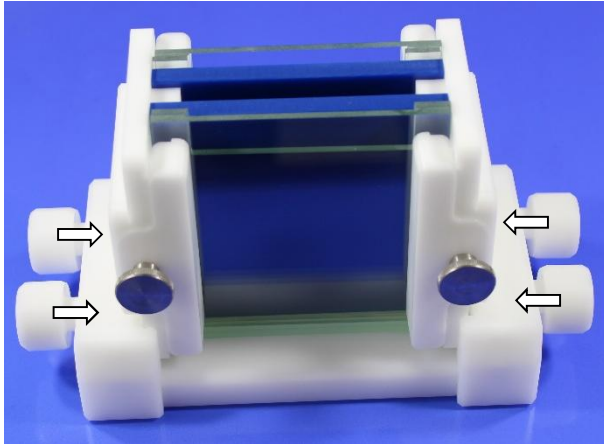
Place the inner core on a flat bench and unloose the screws. Slide in a prepared glass plate sandwich with shorter plate inwards. Take care that the plate edges are in contact with the bench and the sandwich is not wedged. By tightening the screws, the

sandwich is fixed. Turn the inner core and install the second sandwich as described above.

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
Pull the screws of the base unit outwards.

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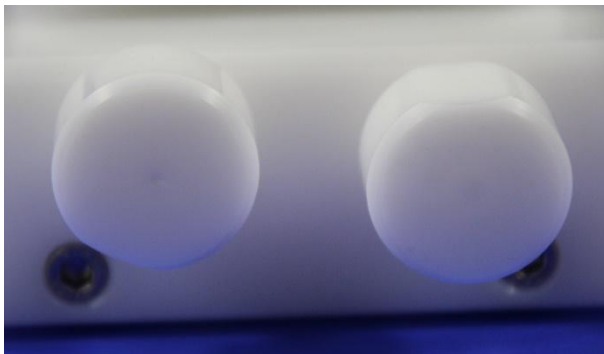
Place the inner core into the base unit and push the screws back inwards.

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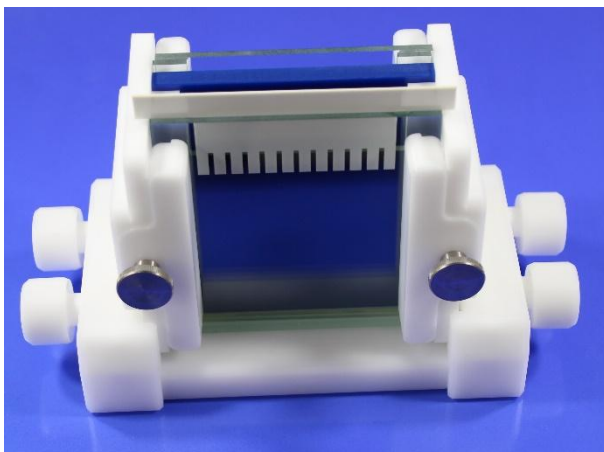Turn the screws crosswise to press the inner core onto the rubber gasket. For better orientation, the screws have a flat side.

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Cast the separation gel. Directly after filling the solution for the stacking gel in, place a comb between the two plates.

**SERVA**  
Electrophoresis

---

**SERVA Electrophoresis GmbH**

**Address** Carl-Benz-Str. 7, 69115 Heidelberg, Germany **Web site** [www.serva.de](http://www.serva.de) **Email** [info@serva.de](mailto:info@serva.de)  
**Customer care Germany** Phone: 06221 13840-46, Fax: 06221 13840-10 **Customer care International**  
Phone: +49 6221 13840-47, Fax: +49 6221 13840-10 **Technical service** Phone: +49 6221 13840-44  
**Technical service Collagenase** Phone: +49 4122 712-413