

Product of the Month

SERVA
■ serving scientists ■

Get **20%** discount* now!

December

*For all SERVA customers within the EU. Outside Germany and Austria online shop orders only. Can not be combined with other discounts, promotions etc.

DNADecon

Highly Efficient against Nucleic Acid Contaminations

DNADecon is a highly effective, ready-to-use spray solution for removing DNA and RNA contaminants on surfaces, equipment and laboratory equipment.

Features

- Ready-to-use
- Non-alkaline
- Non-carcinogenic



Decontamination of surfaces

- Spray **DNADecon** directly onto the surface and clean with a paper towel.
- Rinse with DNA-free water and dry with a clean paper towel.

Decontamination of laboratory devices and porous surfaces

- Apply **DNADecon** on a paper towel and clean the surface, appliances or equipment thoroughly.
- Rinse with DNA-free water and dry with a clean paper towel.

Decontamination of pipettes

- Follow the manufacturer's instructions for pipette disassembling.
- Immerse the stem of the pipette, without the seals, in **DNADecon** for 1 min
- Rinse with DNA-free water, dry and assemble the pipette.

Ordering Information

DNADecon	Cat. no.	Site	List price	Special price*
Spray bottle	39810.01	250 ml	31.00 €	24.80 €
Refill pack	39810.02	500 ml	42.00 €	33.60 €

* The special prices are valid until December 31st, 2020.

For further information,
please contact:

Dr. Judith Koch
Tel.: +49 6221 13840-44
E-Mail: Tech.Service@serva.de

SERVA Electrophoresis GmbH
Carl-Benz-Str. 7
D-69115 Heidelberg • Germany
www.serva.de